

## How to Get Eight Wipes from a Single Pharma-Wipe® Wiper

(This method will substantially reduce your true cost of cleanroom disposables)

Assume: 1) one wipe per surface of the wiper, 2) you are using a 9" x 9" wiper  
3) All folding of wipers is done in mid-air



**Step 1:** Fold the wiper in half, creating a double-thickness 9" x 4½"




**Step 2:** Fold the wiper in half again, creating a four thickness 4½" x 4½"



**Step 3:** Wipe once from the cleanest to the dirtiest area, turn it over and wipe again. Unfold and refold so that the two contaminated sides are together. The result is a four thickness 4½" x 4½" with two clean outer sides exposed and two contaminated sides touching on the inside of the wiper.



**Step 4:** Wipe once, turn it over and wipe again.

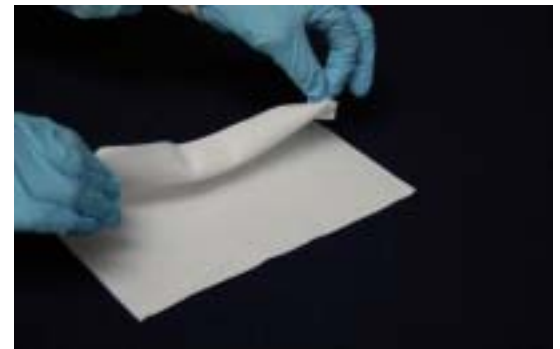

**Step 5:** Unfold the wiper completely and repeat Step 1, ensuring the contaminated sides meet on the initial fold. Repeat Steps 2, 3 & 4.



888-909-7700  
[797@acutecareonline.com](mailto:797@acutecareonline.com)

Pharma-Wipe®