

DISINFECTANTS & CLEANERS



PHARMA-HOL

is a 70% IPA / 30% WFI cleaner that is validated sterile and filtered with a .1 micron filter resulting in extremely low Sub-Visible Particulate levels many times lower than industry specifications.

Dwell time: Leave a surface wet for a minimum of 30 Seconds before wiping with a Pharma-Wipe

PHARMA-D

an EPA registered, one-step disinfectant, cleaner, mildewstat, virucide, fungicide, and sanitizer designed for use in the cleanroom. This USP 797 compliant disinfectant is to be used as a daily disinfectant in rotation with our Pharma-HOL®.

Dwell time: To disinfect leave surface wet for 10 minutes.

DECON-SPORE

an EPA registered Peracetic Acid and Hydrogen Peroxide solution that is for use as a broad-spectrum sanitizer, disinfectant, virucide, fungicide, sporicidal, cleaner and deodorizer in the hospital and healthcare industries.

Dwell Time: Kills Clostridium Difficile Spores in 5 minutes, for all other functions please allow a dwell time of 10 minutes.

For frequency in cleaning see the page below for the cleaning chart.

ROTATION OF CLEANERS & DISINFECTANTS

Site	Cleaning (Pharma-D)	Disinfecting (Pharma-HOL)	Apply Sporidical Disinfectant (Decon-Spore)
PEC(s) and equipment inside the PEC(s)	<ul style="list-style-type: none"> Equipment and all interior surfaces of the PEC daily on days when compounding occurs and when surface contamination is known or suspected 	<ul style="list-style-type: none"> Equipment and all interior surfaces of the PEC daily on days when compounding occurs and when surface contamination is known or suspected 	<ul style="list-style-type: none"> Monthly for entities compounding Category 1 and/or Category 2 CSPs Weekly for entities compounding Category 3 CSPs
Removable work tray of the PEC, when applicable	<ul style="list-style-type: none"> Work surface of the tray daily on days when compounding occurs All surfaces and the area underneath the work tray monthly 	<ul style="list-style-type: none"> Work surface of the tray on days when compounding occurs All surfaces and the area underneath the work tray monthly 	<ul style="list-style-type: none"> Work surfaces of the tray monthly All surfaces and the area underneath the work tray monthly
Pass-Through Chambers	<ul style="list-style-type: none"> Daily on days when compounding occurs 	<ul style="list-style-type: none"> Daily on days when compounding occurs 	<ul style="list-style-type: none"> Monthly for entities compounding Category 1 and/or Category 2 CSPs
Work surface(s) outside the PEC	<ul style="list-style-type: none"> Daily on days when compounding occurs 	<ul style="list-style-type: none"> Daily on days when compounding occurs 	<ul style="list-style-type: none"> Weekly for entities compounding Category 3 CSPs
Floor(s)	<ul style="list-style-type: none"> Daily on days when compounding occurs 	<ul style="list-style-type: none"> Daily on days when compounding occurs 	
Wall(s), door(s), and door frame(s)	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly
Ceiling(s)	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly
Storage shelving and bins	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly
Equipment outside the PEC(s)	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly 	<ul style="list-style-type: none"> Monthly

STANDARDS OF PRACTICE

Cleaning with Pharma-D & Disinfecting with Pharma-HOL in the PEC

- If necessary, remove visible particles, debris, or residue with an appropriate solution (e.g., Sterile Water for Injection or Sterile Water for Irrigation) using a sterile, Pharma-Wipe.
- Using a Sterile Pharma-Wipe, apply Pharma-D and let sit for 10 minutes followed by Pharma-HOL let sit for 30 seconds.
- Using a sterile Pharma-Wipe, apply Pharma-HOL to equipment and all interior surfaces in the PEC.
- Allow the surface to dry completely before beginning compounding.

Procedures for Applying Decon-Spore in the PEC

- If necessary, remove visible particles, debris, or residue with an appropriate solution (e.g., Sterile Water for Injection or Sterile Water for Irrigation) using sterile, Pharma-Wipes.
- After cleaning and disinfecting (see box above), apply the Decon-Spore using a sterile Pharma-Wipe to all surfaces and the area underneath the work tray; if the sporicidal disinfectant is a sterile EPA-registered (or equivalent for entities outside the US) one-step disinfectant sporicidal cleaner, separate cleaning and disinfecting steps are not required.
- Let the applied Decon-Spore sit for 5 minutes.
- Using a sterile Pharma-Wipe, apply sterile 70% IPA to all interior surfaces, including underneath the work tray.
- Allow the surface to dry completely before beginning compounding.

QUESTIONS?
GIVE US A CALL

AcuteCare
PHARMACEUTICALS

A hospenco BRANDS GROUP COMPANY

1-888-909-7700